## UNPACKING


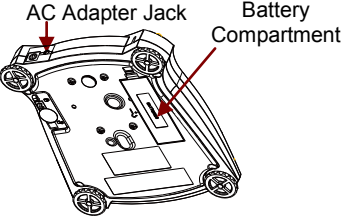
Carefully remove your Scout balance and each of its components from the package. Store the packaging material to ensure safe storage and future transport. Included Components:

- Balance
- Quick Start Guide
- Warranty Card
- AC Adapter & Plug
- Pan
- Stacking & Storage Cover (not applicable for 1mg models)
- Calibration Mass (only for models up to 620g capacity)

## SELECTING THE LOCATION

Place on a firm, steady surface with sufficient space. Avoid excessive air current, vibrations, heat sources, or rapid temperature changes.

## PREPARATION FOR USE
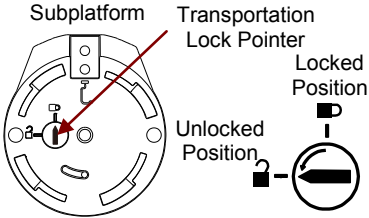
<p><b>Step-1</b></p>	<p><b>Release the red Transportation Lock on the subplatform of the balance</b></p> <p>Turn the pointer 90° counter-clockwise.</p>	
<p><b>Step-2</b></p>	<p><b>Install the Pan</b></p> <p><i>Rectangular pan:</i> Place diagonally on sub-platform and rotate counter-clockwise until it locks.</p> <p><i>Round pan:</i> Set straight down on subplatform.</p>	
<p><b>Step-3</b></p>	<p><b>Connect Power</b></p> <p>Connect the AC Adapter to the balance input jack.</p> <p><i>Battery Installation</i></p> <p>Install the Four "AA" (LR6) batteries (not included) with polarity as shown in the battery compartment.</p>	
<p><b>Step-4</b></p>	<p><b>Level the balance</b></p> <p>Level the balance before using by adjusting the feet until level bubble is center</p>	

## SAFETY PRECAUTIONS

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

## CONTROLS

- **Button Functions**

Zero 	<p><u>Short Press:</u> Turns the balance on</p> <p>If balance is On, sets Zero</p> <p><u>Long Press:</u> Turns the balance Off</p>
Tare	Tare

- **Main Application Screen**

This balance utilizes a touch-sensitive display. *Touch* areas and Buttons to control the balance functions.

<p>Application</p> <p>Stability (*), Net (NET), Gross (G) and/or center of zero (&gt;0&lt;) indicators</p> <p>Reference Fields</p>		<p>Capacity and readability</p> <p>Result Field: Information varies by application</p> <p>Touch <b>g</b> to change unit</p> <p>Application Buttons: Functions vary by application</p>
--	--	---

## CALIBRATION



With balance turned off:

<b>Step-1</b>	Clear Pan.	
<b>Step-2</b>	Turn On balance	
<b>Step-3</b>	Press Menu	
<b>Step-4</b>	Press Calibration	
<b>Step-5</b>	Choose preferred calibration method.	
<b>Step-6</b>	Follow the screen instructions.	

## OPERATION & APPLICATIONS

The balance can be configured to operate in various Application modes. Touch the top left button in the Application area to choose the mode.

The Scout STX Balance incorporates the following Applications:

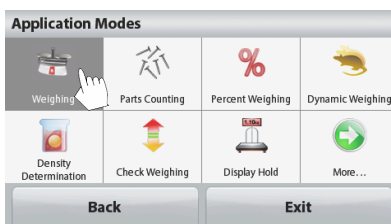
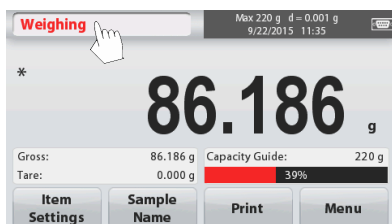
### Weighing

1. Select Weighing (this application is the default).
2. Press **Tare** or **Zero** if necessary to begin.
3. Place objects on the pan to display the weight. When stable, the \* appears. The resulting value is displayed in the main Weighing Line in the active unit of measure.

## MENU & SCREEN NAVIGATION

All menu navigation is performed by touching the display. To enter the Menu, touch **Menu** from any Application Home screen. The Main menu appears, with buttons for **Back** and **Exit**. Continue touching the appropriate list item to navigate to the Menu items or touch a button to change location.

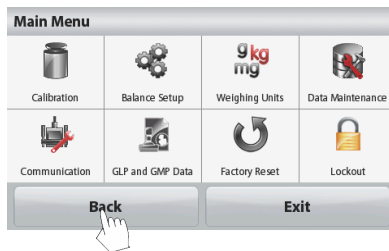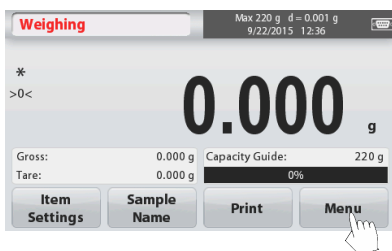**Calibration:**  
Touch to view calibration options.

**Balance Setup:**  
Touch to view balance settings.

**Weighing Units:**  
Touch to view weighing units.

**Data Maintenance:**  
Touch to view data maintenance options.

**Communication:**  
Touch to view COM Device Settings and Print Settings.

**GLP and GMP Data:**  
Insert user data for traceability.

**Factory Reset:**  
Touch to do a Factory reset.

**Lockout:**  
Touch to access the Lockout menu list.


## ADDITIONAL FEATURES

### • **Connecting the Interface**

Use an optional interface connectivity kit to connect the balance either to a computer, printer or OHAUS auxiliary display.

Interface connection on the rear of the balance:



Below Interface kit accessories are available:  
RS232, USB Host, USB Device, Ethernet, Bluetooth\*.

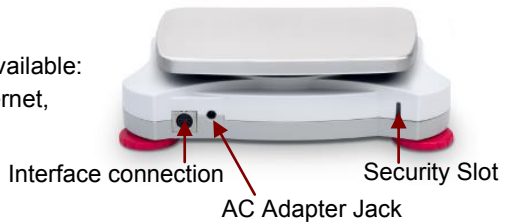\* Interface kits may vary according to local regulations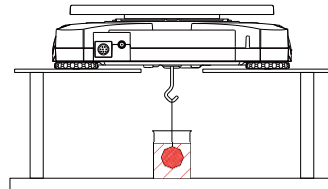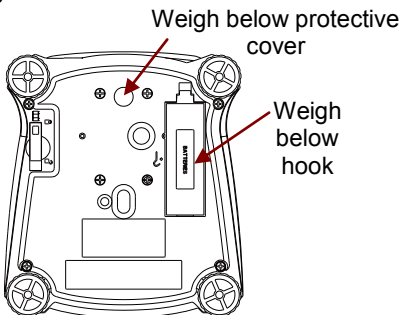

### • **Weigh Below Hook**

The weigh below hook is located at the reverse side of the battery cover as shown. To use this feature, remove the red protective cover underneath for the weigh below opening.

## COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> .
	AS/NZS 61000.6.1, AS/NZS 61000.6.3
	CAN/CSA-C22.2 No. 61010-1 UL Std. No. 61010-1

## INSTRUCTION MANUAL

For more detailed information about the balance, please refer to the product instruction manual. The instruction manual can be obtained either by scanning the QR code or by going to the product page and from there download it.

Go to [www.ohaus.com](http://www.ohaus.com) to register your product's warranty.

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